




# Dilution Bottle Handling Instructions

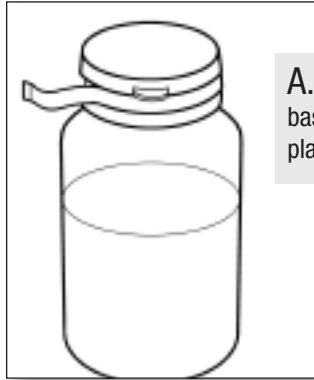

A. Select proper dilution bottle based on cap color. Remove plastic tear-away sterility strip.

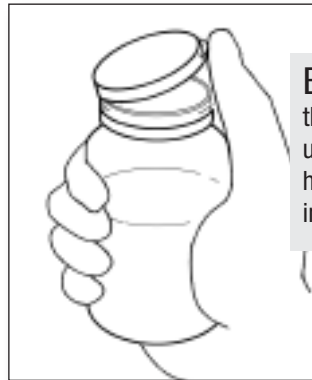


B. To open, grasp bottle in one hand, position thumb directly under cap tab, and carefully push up. Cap will stay in position, allowing use of two hands to add sample to diluent. Do not touch inside of cap or bottle.

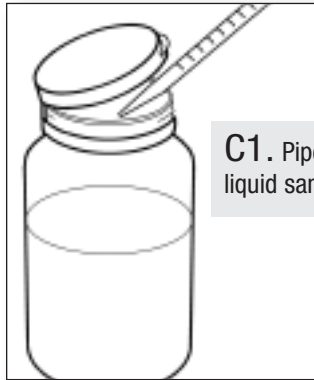

C1. Pipet in liquid samples.

OR

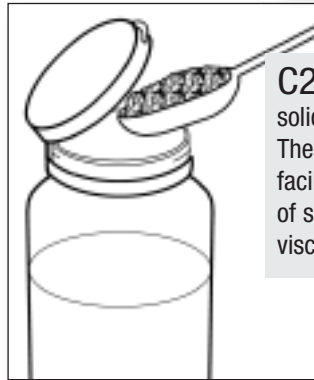

C2. Weigh in solid samples. The wide mouth facilitates addition of solid, bulky, or viscous samples.



D. To close, firmly press down on cap from either behind, using forefingers, or from the front. *Be certain the cap clicks shut all the way around.* To process sample, shake and tap cap once or twice before opening.